

*Bylee*

**Worklist: 1805**

<u>LAB CASE</u>	<u>ITEM</u>	<u>TASK ID</u>	<u>DESCRIPTION</u>	
C2017-1238	1	89728	AM 27 Blood THC Quant by LC	
C2017-1263	1	89729	AM 27 Blood THC Quant by LC	
C2017-1309	1	89730	AM 27 Blood THC Quant by LC	
C2017-1313	1	89731	AM 27 Blood THC Quant by LC	
M2017-2821	2	89732	AM 27 Blood THC Quant by LC	
M2017-3046	2	89733	AM 27 Blood THC Quant by LC	

*A*

# Quantitation of THC and Metabolites in Blood by LC-MS/MS

Extraction Date: 7-11-17

Analyst: Anne Nord

## PRE-ANALYTIC

Plate Lot# Custom - 0490364 Plate Exp. 9-21-2017 External QC Lot 91317, exp 9-13-17

1. Ensure all solutions are within expiration date.
- Mobile Phase A: **0.1% Formic Acid in LCMS Water**
  - Mobile Phase B: **0.1% Formic Acid in LCMS Acetonitrile**
  - **LCMS Methanol**
  - **Blank/Negative Blood: Lot 321632-1**
  - **0.1% Formic Acid in water**
  - **MTBE**
  - **Hexane**
- Column: **UCT Selectra DA 100 x 2.1 mm 3um**
2. Check levels of mobile phases and needle wash and refill as necessary. Ensure waste is not full.
3. Purge Pump and Load appropriate Acq. Method, allow system to equilibrate for approx. 30 min.
4. Create worklist. Data path name: 71117 THC Quant

## ANALYTIC

1. Remove standards plate, blood, and samples from cold storage. Allow to reach room temperature.
2. Add **1000 µL blood** to wells of analytical (standards) plate. Place cover on Plate
- Blank blood for locations containing standards/QCs and internal standards  
Sample blood for locations containing only internal standards
3. Place on shaking incubator at ambient temp., **900rpm for 15 minutes**. *Shaker ID 66759*
4. Pipette **500µL 0.1% formic acid** to all wells of standards plate.
5. Place on shaking incubator at ambient temp., **900rpm for 15 minutes**.
6. Transfer **800µL of blood+acid mixture** to corresponding wells of SLE+ plate.
7. Apply positive pressure for approx. 10-15 seconds (or until no liquid remains on top of sorbent). Wait 5 min. (**Load blood samples at 85- 100 PSI- Selector to Right**)
8. Add **2.25mL MTBE** and allow to flow under gravity for 5 minutes. (**add in 3 increments of 750uL**)
9. Apply positive pressure for approx. 15 seconds (**10-15 PSI- Selector to left -**).
10. Add **2.25mL Hexane** and allow to flow under gravity for 5 minutes. (**add in 3 increments of 750uL**)
11. Apply positive pressure for approx. 15 seconds. (**10-15 PSI Selector to the left**)
12. Remove collection plate containing eluate.
13. Place collection plate on SPE Dry and evaporate to dryness at approx. 35°C. *SPE Dry ID 66819*
14. Reconstitute in **100 µL MeOH** and heat seal plate with foil. Place in autosampler and run worklist.

## POST-ANALYTIC

1. Open quantitation software and create a new quantitation batch.  
Batch name: 71117 cann quant
2. Make any necessary integration changes. Limit curves based on validated linear ranges (3-50ng/mL).
3. Were all appropriate standards used in the curve for each analyte?  Y /  N  
Are  $r^2$  values  $\geq 0.98$  for each analyte?  Y /  N
4. Did all QCs pass for each analyte?  Y /  N Were QCs entered into QC charting?  Y /  N
5. Central File Packet to include:  LIMS Worklist:  Method Checklist  Calibration and Control Reports

COMMENTS

# ISP FORENSICS - Cd'A Instrument # 62340


## Cannabinoids Analysis Report

**Batch Data Path** D:\2017 Data\71117 THC Quant\QuantResults\71117 cann quant.batch.bin  
**Analysis Time** 7/12/2017 3:38 PM **Analyst Name** ISP Tox  
**Report Time** 7/12/2017 3:40 PM **Reporter Name** ISP Tox  
**Last Calib Update** 7/12/2017 3:38 PM **Batch State** Processed

### Analysis Info

**Acq Time** 2017-07-12 12:12 **Data File** Neg Control.d  
**Sample Type** Sample **Sample Name** Neg Control  
**Dilution** 1 **Acq Method** Quant THC 2017.m  
**Position** P2-a2 **Sample Info**  
**Inj Vol** -1 **Comment** AM 27 cannabinoid confirmation

### Sample Chromatogram

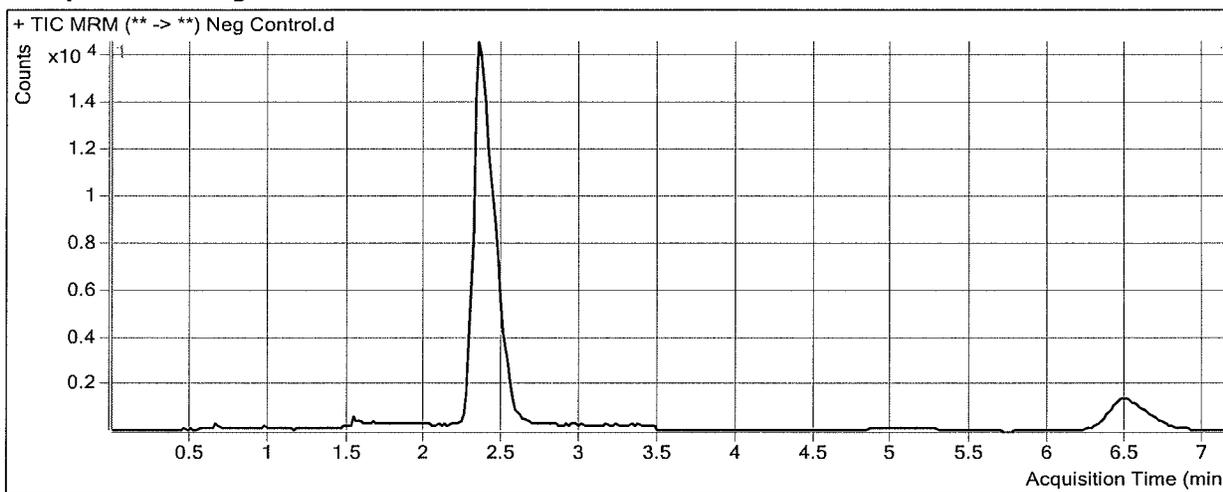

# ISP FORENSICS - Cd'A Instrument # 62340


## Cannabinoids Analysis Report

**Batch Data Path** D:\2017 Data\71117 THC Quant\QuantResults\71117 cann quant.batch.bin  
**Analysis Time** 7/12/2017 3:38 PM **Analyst Name** ISP Tox  
**Report Time** 7/12/2017 3:40 PM **Reporter Name** ISP Tox  
**Last Calib Update** 7/12/2017 3:38 PM **Batch State** Processed

### Analysis Info

**Acq Time** 2017-07-12 12:24 **Data File** QC 10.d  
**Sample Type** QC **Sample Name** QC 10  
**Dilution** 1 **Acq Method** Quant THC 2017.m  
**Position** P2-H1 **Sample Info**  
**Inj Vol** -1 **Comment** AM 27 cannabinoid confirmation

### Sample Chromatogram

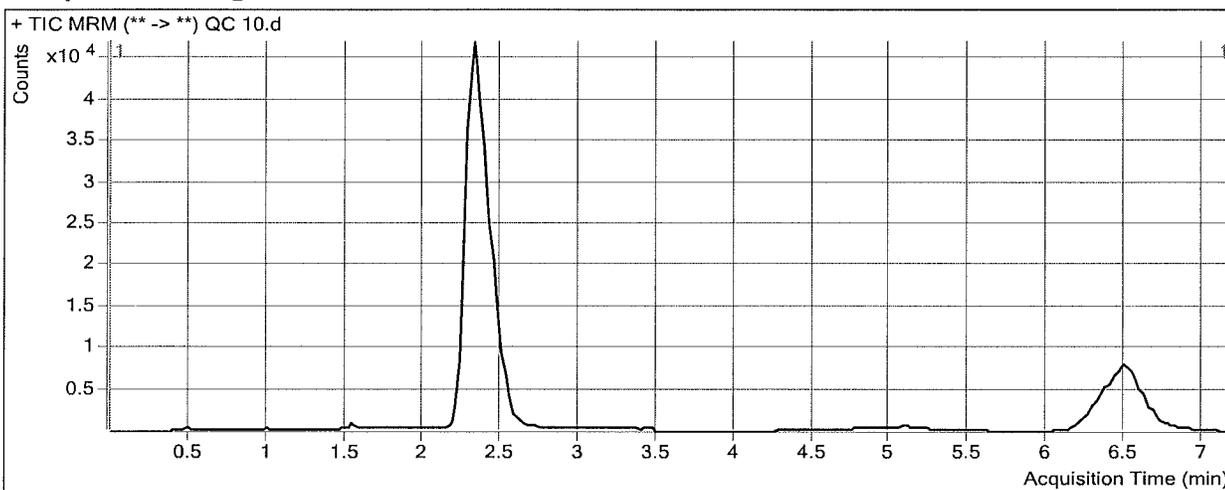

### Results

Compound	ISTD Compound	RT	Response	ISTD Resp	Resp Ratio	Final Conc
THC-OH	THC-OH-d3	2.336	28585	342610	0.0834	9.6590
THC-COOH	THC-COOH-d9	2.486	20666	104303	0.1981	10.0967
THC	THC-d3	6.493	16165	132360	0.1221	10.4661

# ISP FORENSICS - Cd'A Instrument # 62340


## Cannabinoids Analysis Report

**Batch Data Path** D:\2017 Data\71117 THC Quant\QuantResults\71117 cann quant.batch.bin  
**Analysis Time** 7/12/2017 3:38 PM **Analyst Name** ISP Tox  
**Report Time** 7/12/2017 3:40 PM **Reporter Name** ISP Tox  
**Last Calib Update** 7/12/2017 3:38 PM **Batch State** Processed

### Analysis Info

**Acq Time** 2017-07-12 12:36 **Data File** External control lot 91317.d  
**Sample Type** Sample **Sample Name** External control lot 91317  
**Dilution** 1 **Acq Method** Quant THC 2017.m  
**Position** p2b2 **Sample Info**  
**Inj Vol** -1 **Comment** AM 27 cannabinoid confirmation 10 ng

### Sample Chromatogram

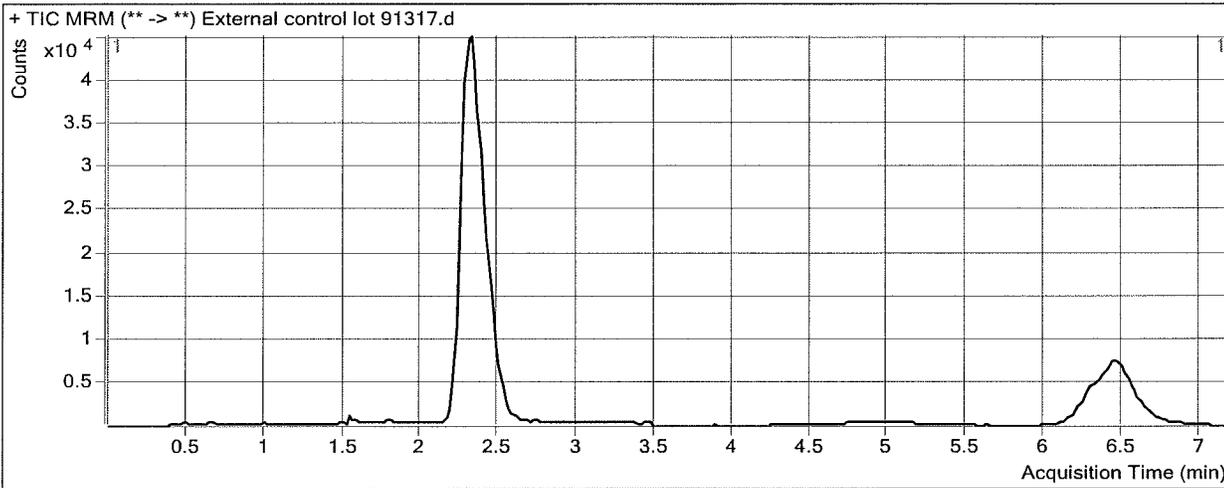

### Results

Compound	ISTD Compound	RT	Response	ISTD Resp	Resp Ratio	Final Conc
THC-OH	THC-OH-d3	2.336	29795	337049	0.0884	10.2285
THC-COOH	THC-COOH-d9	2.466	16324	102008	0.1600	7.9902
THC	THC-d3	6.433	11786	103655	0.1137	9.7437

# ISP FORENSICS - Cd'A Instrument # 62340


## Cannabinoids Analysis Report

**Batch Data Path** D:\2017 Data\71117 THC Quant\QuantResults\71117 cann quant.batch.bin  
**Analysis Time** 7/12/2017 3:38 PM **Analyst Name** ISP Tox  
**Report Time** 7/12/2017 3:40 PM **Reporter Name** ISP Tox  
**Last Calib Update** 7/12/2017 3:38 PM **Batch State** Processed

### Analysis Info

**Acq Time** 2017-07-12 10:37 **Data File** Cal 1- 3ng.d  
**Sample Type** Calibration **Sample Name** Cal 1- 3ng  
**Dilution** 1 **Acq Method** Quant THC 2017.m  
**Position** P2-A1 **Sample Info**  
**Inj Vol** -1 **Comment** AM 27 cannabinoid confirmation

### Sample Chromatogram

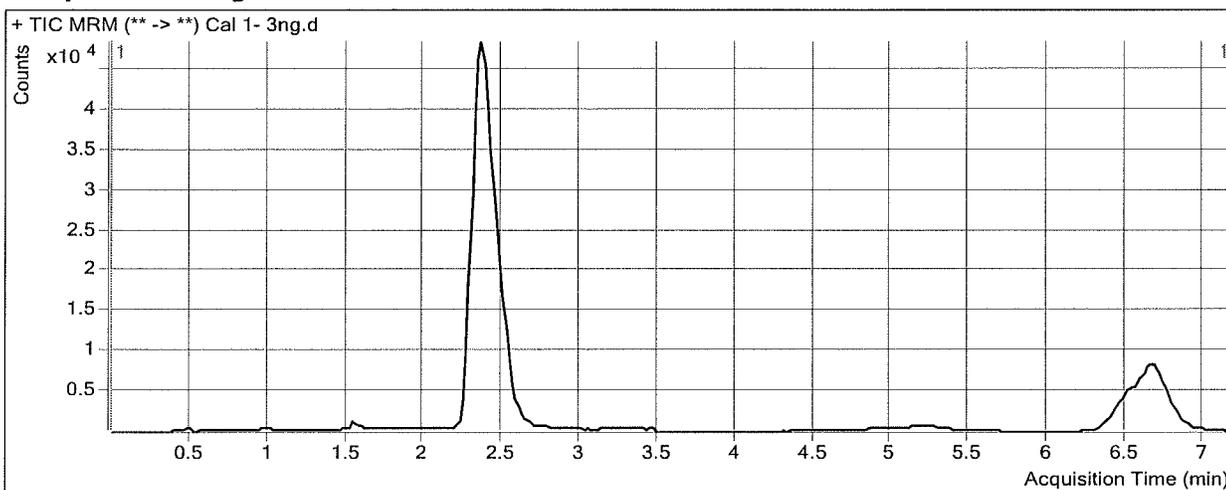

### Results

Compound	ISTD Compound	RT	Response	ISTD Resp	Resp Ratio	Final Conc
THC-OH	THC-OH-d3	2.376	9682	385914	0.0251	2.9716
THC-COOH	THC-COOH-d9	2.526	7975	116503	0.0685	2.9281
THC	THC-d3	6.693	4834	149246	0.0324	2.7696

# ISP FORENSICS - Cd'A Instrument # 62340


## Cannabinoids Analysis Report

**Batch Data Path** D:\2017 Data\71117 THC Quant\QuantResults\71117 cann quant.batch.bin  
**Analysis Time** 7/12/2017 3:38 PM **Analyst Name** ISP Tox  
**Report Time** 7/12/2017 3:40 PM **Reporter Name** ISP Tox  
**Last Calib Update** 7/12/2017 3:38 PM **Batch State** Processed

### Analysis Info

**Acq Time** 2017-07-12 10:49 **Data File** cal 2- 5ng.d  
**Sample Type** Calibration **Sample Name** cal 2- 5ng  
**Dilution** 1 **Acq Method** Quant THC 2017.m  
**Position** P2-B1 **Sample Info**  
**Inj Vol** -1 **Comment** AM 27 cannabinoid confirmation

### Sample Chromatogram

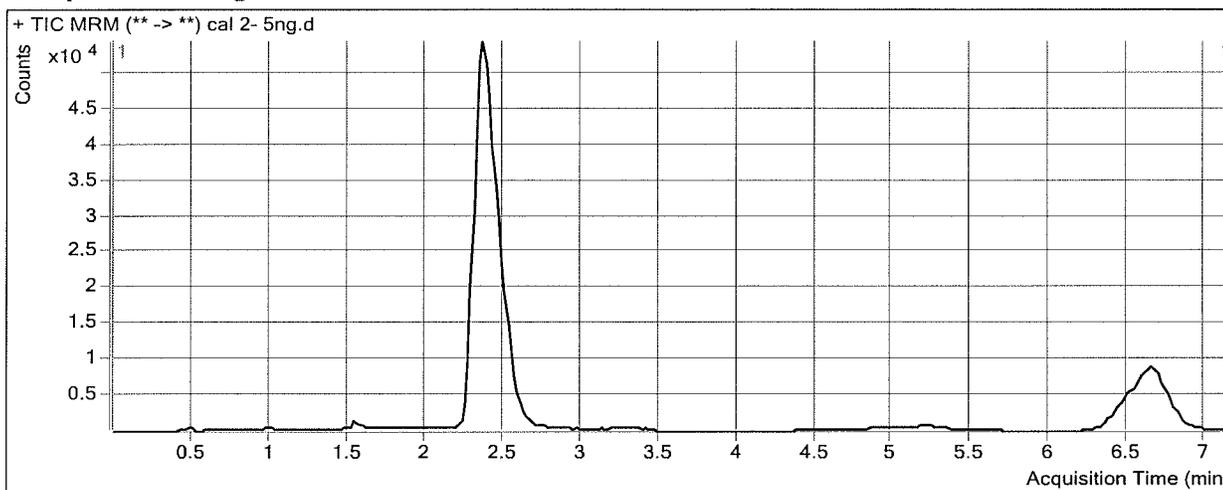

### Results

Compound	ISTD Compound	RT	Response	ISTD Resp	Resp Ratio	Final Conc
THC-OH	THC-OH-d3	2.396	17973	432274	0.0416	4.8616
THC-COOH	THC-COOH-d9	2.526	12906	123475	0.1045	4.9220
THC	THC-d3	6.653	9166	154997	0.0591	5.0633

# ISP FORENSICS - Cd'A Instrument # 62340


## Cannabinoids Analysis Report

**Batch Data Path** D:\2017 Data\71117 THC Quant\QuantResults\71117 cann quant.batch.bin  
**Analysis Time** 7/12/2017 3:38 PM **Analyst Name** ISP Tox  
**Report Time** 7/12/2017 3:40 PM **Reporter Name** ISP Tox  
**Last Calib Update** 7/12/2017 3:38 PM **Batch State** Processed

### Analysis Info

**Acq Time** 2017-07-12 11:01 **Data File** cal 3 - 10ng.d  
**Sample Type** Calibration **Sample Name** cal 3 - 10ng  
**Dilution** 1 **Acq Method** Quant THC 2017.m  
**Position** P2-C1 **Sample Info**  
**Inj Vol** -1 **Comment** AM 27 cannabinoid confirmation

### Sample Chromatogram

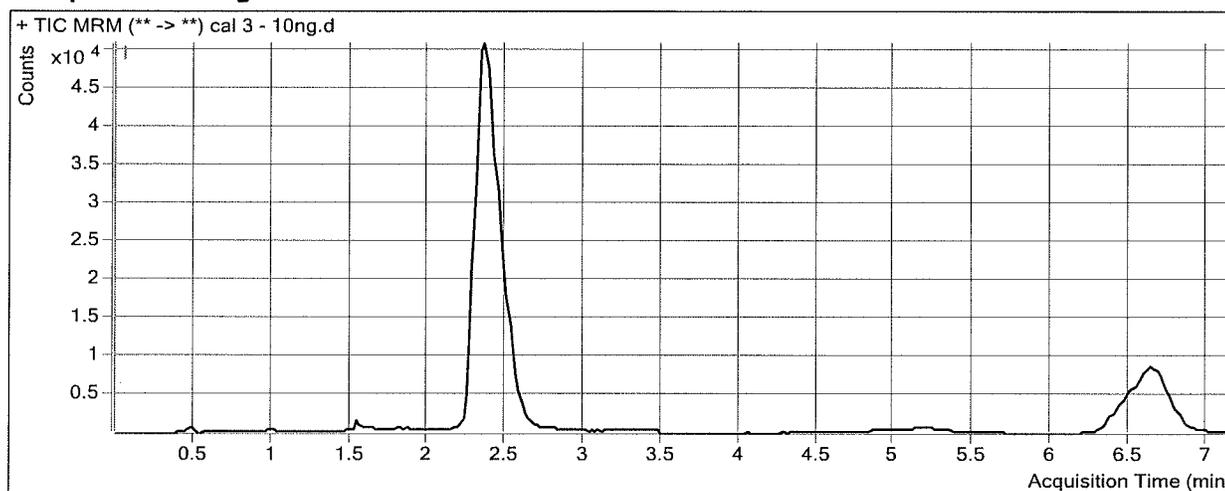

### Results

Compound	ISTD Compound	RT	Response	ISTD Resp	Resp Ratio	Final Conc
THC-OH	THC-OH-d3	2.376	32754	385051	0.0851	9.8462
THC-COOH	THC-COOH-d9	2.506	22399	113451	0.1974	10.0579
THC	THC-d3	6.633	16861	143263	0.1177	10.0855

# ISP FORENSICS - Cd'A Instrument # 62340


## Cannabinoids Analysis Report

**Batch Data Path** D:\2017 Data\71117 THC Quant\QuantResults\71117 cann quant.batch.bin  
**Analysis Time** 7/12/2017 3:38 PM **Analyst Name** ISP Tox  
**Report Time** 7/12/2017 3:40 PM **Reporter Name** ISP Tox  
**Last Calib Update** 7/12/2017 3:38 PM **Batch State** Processed

### Analysis Info

**Acq Time** 2017-07-12 11:13 **Data File** cal 4 - 25ng.d  
**Sample Type** Calibration **Sample Name** cal 4 - 25ng  
**Dilution** 1 **Acq Method** Quant THC 2017.m  
**Position** P2-D1 **Sample Info**  
**Inj Vol** -1 **Comment** AM 27 cannabinoid confirmation

### Sample Chromatogram

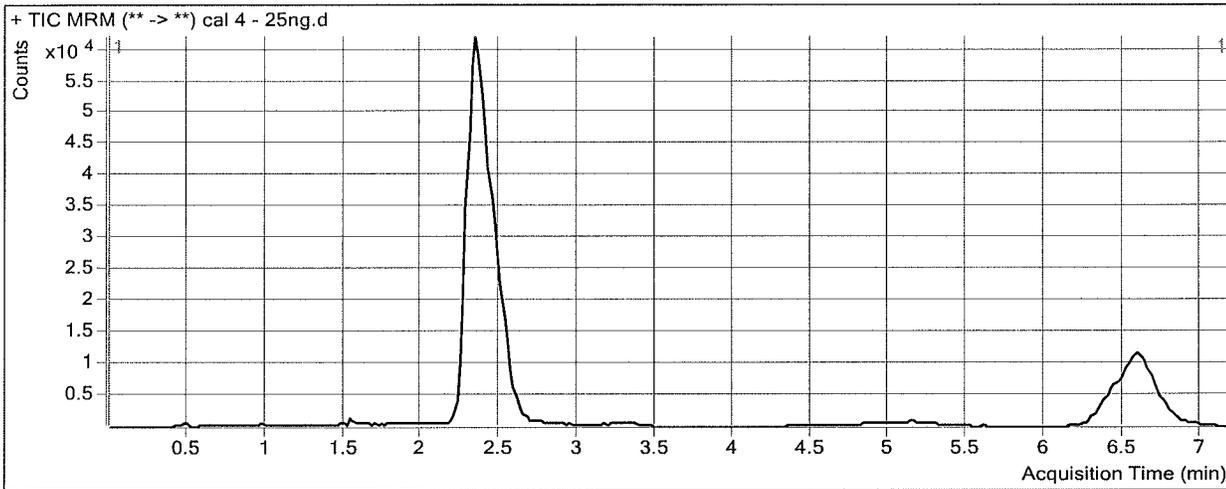

### Results

Compound	ISTD Compound	RT	Response	ISTD Resp	Resp Ratio	Final Conc
THC-OH	THC-OH-d3	2.356	88638	412568	0.2148	24.7218
THC-COOH	THC-COOH-d9	2.506	57270	124327	0.4606	24.6076
THC	THC-d3	6.593	46882	159796	0.2934	25.1549

# ISP FORENSICS - Cd'A Instrument # 62340


## Cannabinoids Analysis Report

**Batch Data Path** D:\2017 Data\71117 THC Quant\QuantResults\71117 cann quant.batch.bin  
**Analysis Time** 7/12/2017 3:38 PM **Analyst Name** ISP Tox  
**Report Time** 7/12/2017 3:40 PM **Reporter Name** ISP Tox  
**Last Calib Update** 7/12/2017 3:38 PM **Batch State** Processed

### Analysis Info

**Acq Time** 2017-07-12 11:25 **Data File** cal 5 - 50ng.d  
**Sample Type** Calibration **Sample Name** cal 5 - 50ng  
**Dilution** 1 **Acq Method** Quant THC 2017.m  
**Position** P2-E1 **Sample Info**  
**Inj Vol** -1 **Comment** AM 27 cannabinoid confirmation

### Sample Chromatogram

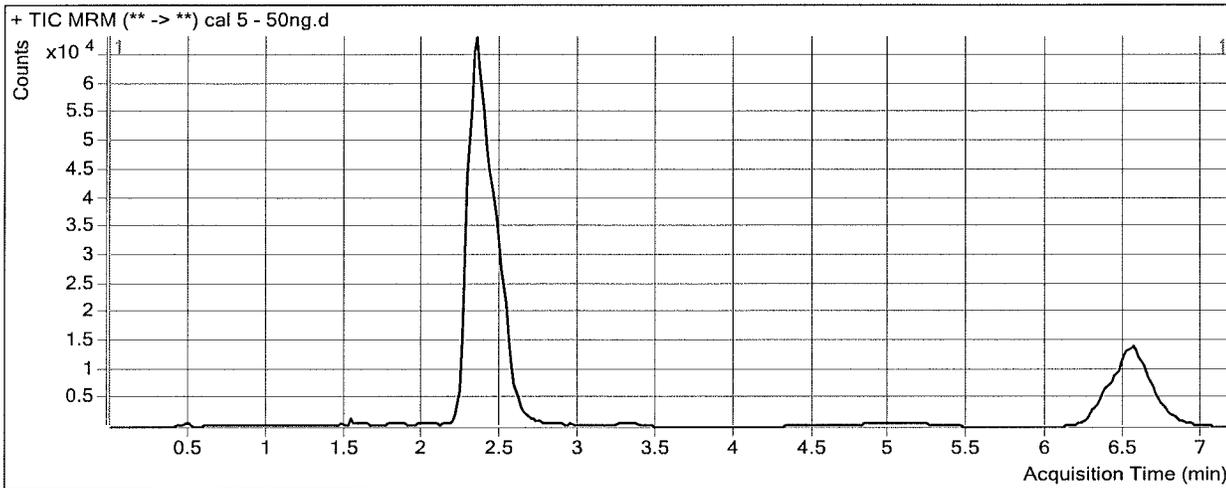

### Results

Compound	ISTD Compound	RT	Response	ISTD Resp	Resp Ratio	Final Conc
THC-OH	THC-OH-d3	2.356	175095	364386	0.4805	55.1743
THC-COOH	THC-COOH-d9	2.486	111217	109721	1.0136	55.1760
THC	THC-d3	6.553	89864	140601	0.6391	54.8099

# ISP FORENSICS - Cd'A Instrument # 62340


## Cannabinoids Analysis Report

**Batch Data Path** D:\2017 Data\71117 THC Quant\QuantResults\71117 cann quant.batch.bin  
**Analysis Time** 7/12/2017 3:38 PM **Analyst Name** ISP Tox  
**Report Time** 7/12/2017 3:40 PM **Reporter Name** ISP Tox  
**Last Calib Update** 7/12/2017 3:38 PM **Batch State** Processed

### Analysis Info

**Acq Time** 2017-07-12 11:37 **Data File** cal 6 - 100ng.d  
**Sample Type** Calibration **Sample Name** cal 6 - 100ng  
**Dilution** 1 **Acq Method** Quant THC 2017.m  
**Position** P2-F1 **Sample Info**  
**Inj Vol** -1 **Comment** AM 27 cannabinoid confirmation

### Sample Chromatogram

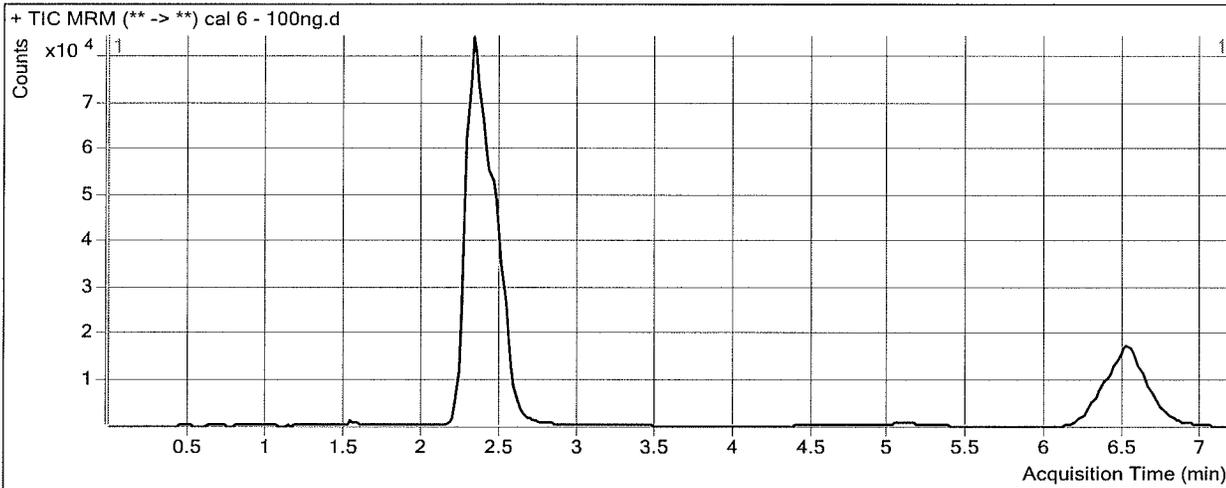

### Results

Compound	ISTD Compound	RT	Response	ISTD Resp	Resp Ratio	Final Conc
THC-OH	THC-OH-d3	2.336	300287	357366	0.8403	96.4102
THC-COOH	THC-COOH-d9	2.466	181663	105732	1.7181	94.1206
THC	THC-d3	6.513	148401	133329	1.1130	95.4556

# ISP FORENSICS - Cd'A Instrument # 62340


## Cannabinoids Analysis Report

**Batch Data Path** D:\2017 Data\71117 THC Quant\QuantResults\71117 cann quant.batch.bin  
**Analysis Time** 7/12/2017 3:38 PM **Analyst Name** ISP Tox  
**Report Time** 7/12/2017 3:40 PM **Reporter Name** ISP Tox  
**Last Calib Update** 7/12/2017 3:38 PM **Batch State** Processed

### Analysis Info

**Acq Time** 2017-07-12 11:48 **Data File** cal 7 - 250ng.d  
**Sample Type** Calibration **Sample Name** cal 7 - 250ng  
**Dilution** 1 **Acq Method** Quant THC 2017.m  
**Position** P2-G1 **Sample Info**  
**Inj Vol** -1 **Comment** AM 27 cannabinoid confirmation

### Sample Chromatogram

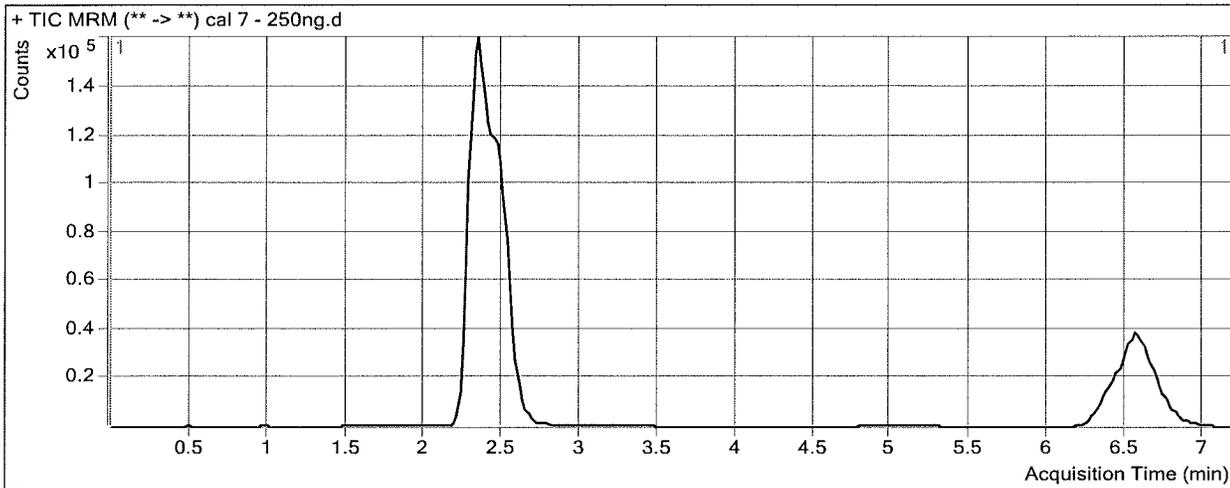

### Results


Compound	ISTD Compound	RT	Response	ISTD Resp	Resp Ratio	Final Conc
THC-OH	THC-OH-d3	2.356	859525	395794	2.1716	249.0144
THC-COOH	THC-COOH-d9	2.486	508829	111597	4.5595	251.1878
THC	THC-d3	6.553	427941	147010	2.9110	249.6612

# ISP Forensics Calibration Curve Report

**Batch Data Path** D:\2017 Data\71117 THC Quant\QuantResults\71117 cann quant.batch.bin

**Last Calib Update** 7/12/2017 3:38 PM **Analyst Name** ISP TOX

**Target Compound** *THC-OH*  
**Internal Standard** *THC-OH-d3*

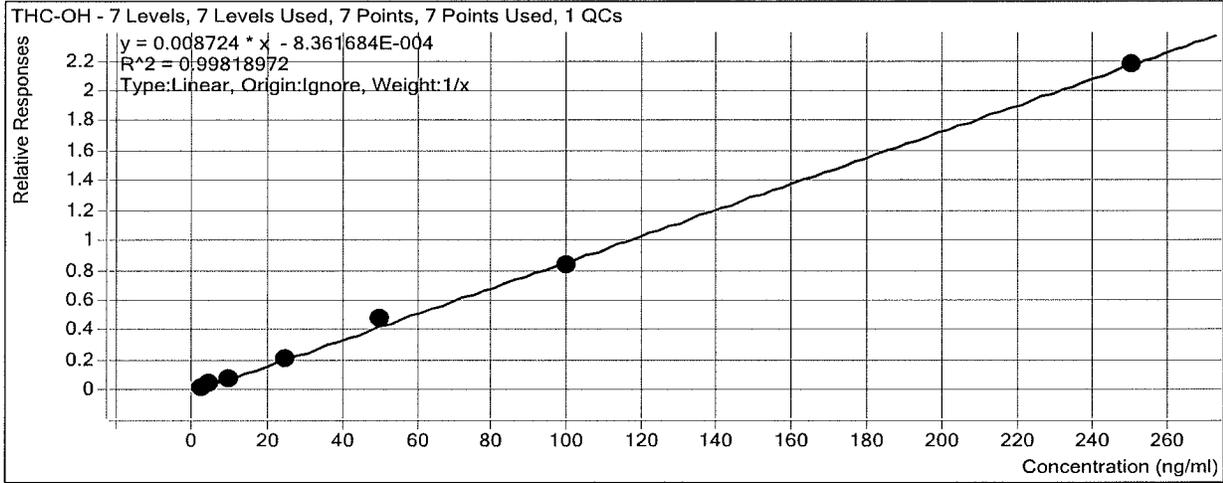


Sample	Level	Enabled	Exp Conc	Final Conc	Accuracy
Cal 1- 3ng	1	<input checked="" type="checkbox"/>	3	3.0	99.1
cal 2- 5ng	2	<input checked="" type="checkbox"/>	5	4.9	97.2
cal 3 - 10ng	3	<input checked="" type="checkbox"/>	10	9.8	98.5
QC 10	3	<input checked="" type="checkbox"/>	10	9.7	96.6
cal 4 - 25ng	4	<input checked="" type="checkbox"/>	25	24.7	98.9
cal 5 - 50ng	5	<input checked="" type="checkbox"/>	50	55.2	110.3
cal 6 - 100ng	6	<input checked="" type="checkbox"/>	100	96.4	96.4
cal 7 - 250ng	7	<input checked="" type="checkbox"/>	250	249.0	99.6

# ISP Forensics Calibration Curve Report

**Batch Data Path** D:\2017 Data\71117 THC Quant\QuantResults\71117 cann quant.batch.bin

**Last Calib Update** 7/12/2017 3:38 PM **Analyst Name** ISP TOX

**Target Compound** *THC-COOH*  
**Internal Standard** *THC-COOH-d9*

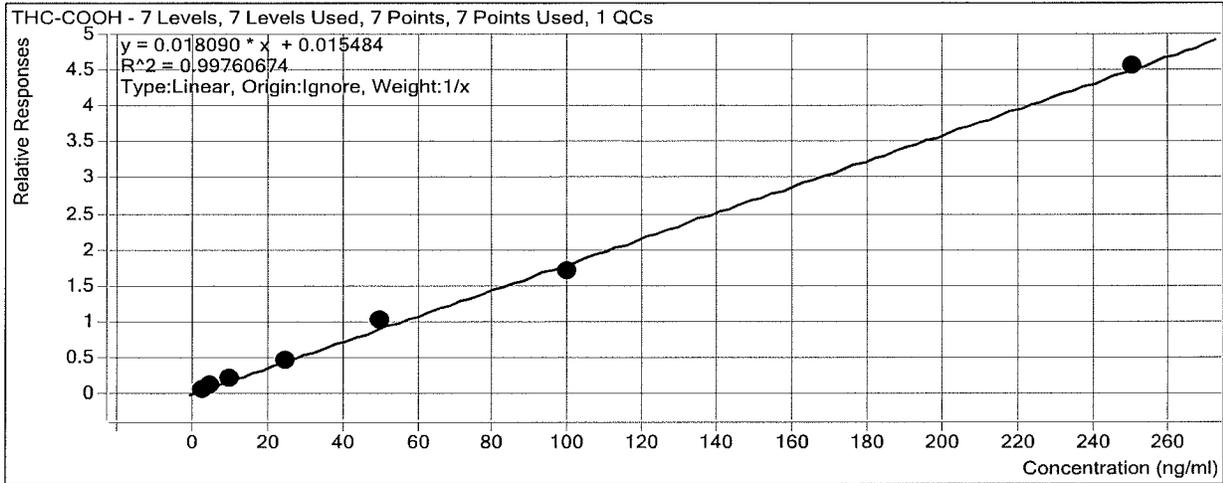


Sample	Level	Enabled	Exp Conc	Final Conc	Accuracy
Cal 1- 3ng	1	<input checked="" type="checkbox"/>	3	2.9	97.6
cal 2- 5ng	2	<input checked="" type="checkbox"/>	5	4.9	98.4
cal 3 - 10ng	3	<input checked="" type="checkbox"/>	10	10.1	100.6
QC 10	3	<input checked="" type="checkbox"/>	10	10.1	101.0
cal 4 - 25ng	4	<input checked="" type="checkbox"/>	25	24.6	98.4
cal 5 - 50ng	5	<input checked="" type="checkbox"/>	50	55.2	110.4
cal 6 - 100ng	6	<input checked="" type="checkbox"/>	100	94.1	94.1
cal 7 - 250ng	7	<input checked="" type="checkbox"/>	250	251.2	100.5

# ISP Forensics Calibration Curve Report

**Batch Data Path**                    D:\2017 Data\71117 THC Quant\QuantResults\71117 cann quant.batch.bin

**Last Calib Update**                7/12/2017 3:38 PM                    **Analyst Name**                    ISP TOX

**Target Compound**                *THC*  
**Internal Standard**                *THC-d3*

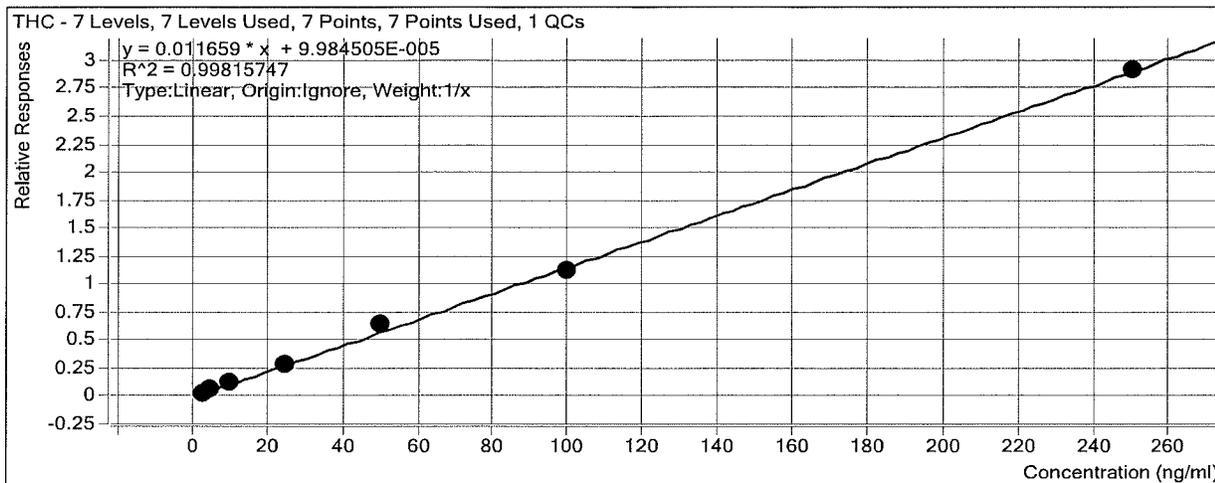

Sample	Level	Enabled	Exp Conc	Final Conc	Accuracy
Cal 1- 3ng	1	<input checked="" type="checkbox"/>	3	2.8	92.3
cal 2- 5ng	2	<input checked="" type="checkbox"/>	5	5.1	101.3
cal 3 - 10ng	3	<input checked="" type="checkbox"/>	10	10.1	100.9
QC 10	3	<input checked="" type="checkbox"/>	10	10.5	104.7
cal 4 - 25ng	4	<input checked="" type="checkbox"/>	25	25.2	100.6
cal 5 - 50ng	5	<input checked="" type="checkbox"/>	50	54.8	109.6
cal 6 - 100ng	6	<input checked="" type="checkbox"/>	100	95.5	95.5
cal 7 - 250ng	7	<input checked="" type="checkbox"/>	250	249.7	99.9